


Specification

AEBSF Hydrochloride *BioChemica*

A1421

Solubility:	50 g/L (H ₂ O)
Physical Description:	Solid
Product Code:	A1421
Product Name:	AEBSF Hydrochloride <i>BioChemica</i>
Specifications:	Assay (titr., calc. on dried substance): min. 98 % Assay (HPLC): min. 98 % Loss on drying: max. 1.0 %
Hazard pictograms	
WGK:	1
Storage:	2-8°C
Signal Word:	Danger
GHS Symbols:	GHS05
H Phrases:	H314
P Phrases:	P280 P305+P351+P338 P310
Molecular Formula:	C ₈ H ₁₀ FNO ₂ S · HCl
M:	239.69 g/mol
CAS:	30827-99-7
CS:	29214900

AppliChem GmbH

Ottoweg 4 • D-64291 Darmstadt • Phone +49 6151 9357 0 • Fax +49 6151 9357 11 • info.de@itwreagents.com • www.itwreagents.com
CEO Joan Roget • Commerzbank Darmstadt • Bank 508 800 50 • Account 0186989900 IBAN DE24 5088 0050 0186 9899 00 • Swiftcode DRESDEFF508 • Finanzamt Darmstadt 07 228 16476 • Register court Darmstadt HRB Nr. 7340

Specification

AEBSF Hydrochloride BioChemica

A1421

Comment

AEBSF irreversibly inactivates thrombin and other serine proteases (e. g. chymotrypsin, kallikrein, plasmin, proteinase K, trypsin) by sulfonylation of a functional group in the active center of the enzyme (1, 2). AEBSF (LD₅₀ oral, mouse 2.8 g/kg) is a substantially less toxic substitute for PMSF and DFP. It is highly water-soluble and its solutions have an acidic pH. At this pH, stability is high. AEBSF is inactivated at 37°C by 50 % after 5 hours. We recommend storage of the stock solution (e. g. 20 mM or 100 mM in water or buffer) at -20°C (stable for up to 2 months; at +4°C approx. 1 week; ref. 6). It should be added to the assay buffers just before use at a final concentration of 0.1 - 2 mM (e. g. ref. 3-5).

Bibliography

(1)Walsmann, P. *et al.* (1972) *Acta biol. med. germ.* **28**, 577-585Inactivation of Trypsin and Thrombin through 4-Amidinobenzene sulfonyl fluoride and 4-(2-Aminoethyl)-benzene sulfonyl fluoride. (2)Markwardt, F. *et al.* (1973) *Thrombosis Res.* **2**, 343-348Synthetic low molecular weight inhibitors of serum kallikrein. (3)Murphy, B.J. *et al.* (1993) *J. Biol. Chem.* **268**, 27355-27362Identification of the sites of selective phosphorylation and dephosphorylation of the rat brain Na⁺ channel a subunit by cAMP-dependent protein kinase and phosphoprotein phosphatases. (4)Nathan, D.F. & Lindquist, S. (1995) *Mol. Cell. Biol.* **15**, 3917-3925Mutational analysis of HSP90 function\: Interactions with a steroid receptor and a protein kinase. (5)Taylor, J.A. *et al.* (1995) *Mol. Cell. Biol.* **15**, 4149-4157Activation of the high-affinity immunoglobulin E receptor FcεRI in RBL-2H3 cells is inhibited by syk SH2 domains. (6)Taylor, J.A. *et al.* (1995) *Immunology* **86**, 629-635Serine protease inhibitors block the priming of monocytes for enhanced superoxide release.

AppliChem GmbH

Ottoweg 4 • D-64291 Darmstadt • Phone +49 6151 9357 0 • Fax +49 6151 9357 11 • info.de@itwreagents.com • www.itwreagents.com
CEO Joan Roget • Commerzbank Darmstadt • Bank 508 800 50 • Account 0186989900 IBAN DE24 5088 0050 0186 9899 00 • Swiftcode DRESDEFF508 • Finanzamt Darmstadt 07 228 16476 • Register court Darmstadt HRB Nr. 7340